

# University of Minnesota Nano Fabrication Center

## Standard Operating Procedure

<b>Equipment name:</b>	Xetch e1 series	<b>Model:</b>	e1
<b>Badger name:</b>	Xef <sub>2</sub> -etch	<b>Revisionist:</b>	Paul Kimani
<b>Location:</b>	Bay 3	<b>Date:</b>	April 25, 2013
<b>Revision #:</b>	1		



### Introduction

This is a system designed to expose samples to xenon di-fluoride (XeF<sub>2</sub>) gas in a cyclic or pulsed mode. XeF<sub>2</sub> can be used to isotropically etch Si, Mo, Ge, Nb, polysilicon, Ti and W. XeF<sub>2</sub> can be used to make very long undercuts with little or no degradation of etch stop, mask or device layers. Advantages of using this system include:

- *Selectivity during etching:* XeF<sub>2</sub> shows very high selectivity vs. silicon to photoresist, SiO<sub>2</sub>, Si<sub>3</sub>N<sub>4</sub>, and aluminum. Selectivity of silicon nitride is better than 100:1 and the selectivity to silicon dioxide is better than 1000:1. No external energy or ion bombardment takes place, and this is the reason XeF<sub>2</sub> exhibits high selectivity to many metals, dielectrics and polymers. The etch reaction is exothermic and results in an increase in chamber temperature. A **short dip in BOE** is useful to break through the oxide.
- *No release stiction:* XeF<sub>2</sub> etching is a dry process so no drying is needed which avoids the sticking issues that often plague wet release processes. Since XeF<sub>2</sub> etching is a dry, room temperature process, delicate structures can be released.
- The system can etch silicon from small part to 100mm (4 inch) wafers. The etch rate is dependent on the surface area of the silicon and decreases with increasing area.

# University of Minnesota Nano Fabrication Center

## Standard Operating Procedure

### Safety

- XeF<sub>2</sub> gas will form HF in presence of water vapor therefore proper ventilation and standard acid precautions are required when handling the bottle exchange (this is a maintenance task)

### Restrictions/Requirements

- Ensure samples are fairly dry of moisture. Bake of sample prior to loading in the etch chamber is recommended to reduce moisture
- Must be qualified to use the tool
- Do not use standard tape with this tool. Kapton or any teflon based tape is okay to use

### Required facilities

- Dry compressed air (70 – 100 psi)
- Vacuum
- Nitrogen gas
- 120V AC Electrical power
- Fume and pump exhaust
- XeF<sub>2</sub> cylinder
- Computer and screen

### Definitions

- Xenon di-fluoride (XeF<sub>2</sub>): a white crystalline substance resembling rock salt. At room temperature, the vapor pressure of XeF<sub>2</sub> is 3.9 torr.
- Expansion chamber: Where sublimated XeF<sub>2</sub> and nitrogen collect before entering the process chamber. Note that N<sub>2</sub> is used for venting, purging and is also combined with XeF<sub>2</sub> for etching.
- Process chamber: Where etching occurs for the amount of time specified by the user.
- Vacuum pump: pumps gases out of both the process and expansion chambers.
- Shower head stop:

### Operation

- If the system has been off, press the ON (green) button to start the e1 series and the vacuum pump. The system has to warm up for four (4) hours before proceeding. If system is on, proceed to (e) below.




- Turn on the computer system (if it has been off)
- The Xetch program will auto start; if it doesn't not Press START- PROGRAMS- XACTIX- Xetch to start it manually.

# University of Minnesota Nano Fabrication Center

## Standard Operating Procedure

- d. After initialization, a dialog box will appear. Press NORMAL startup and allow routine to complete.

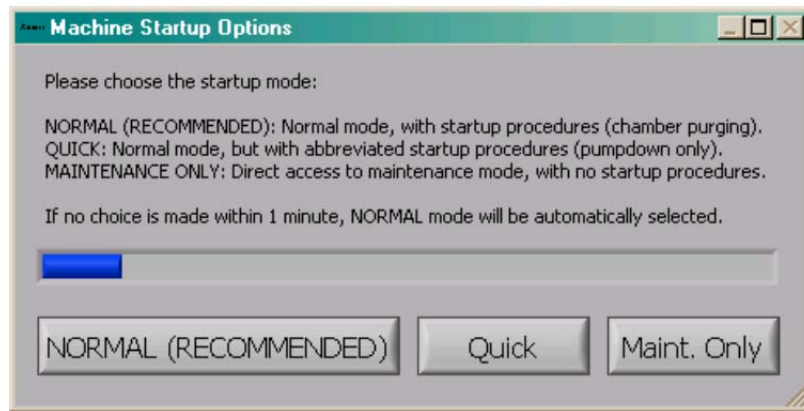

- e. Enter username and password to proceed. Username is USER and password is 1234.




# University of Minnesota Nano Fabrication Center

## Standard Operating Procedure

### Sample loading

- Upon successful login, the main menu below will be displayed.

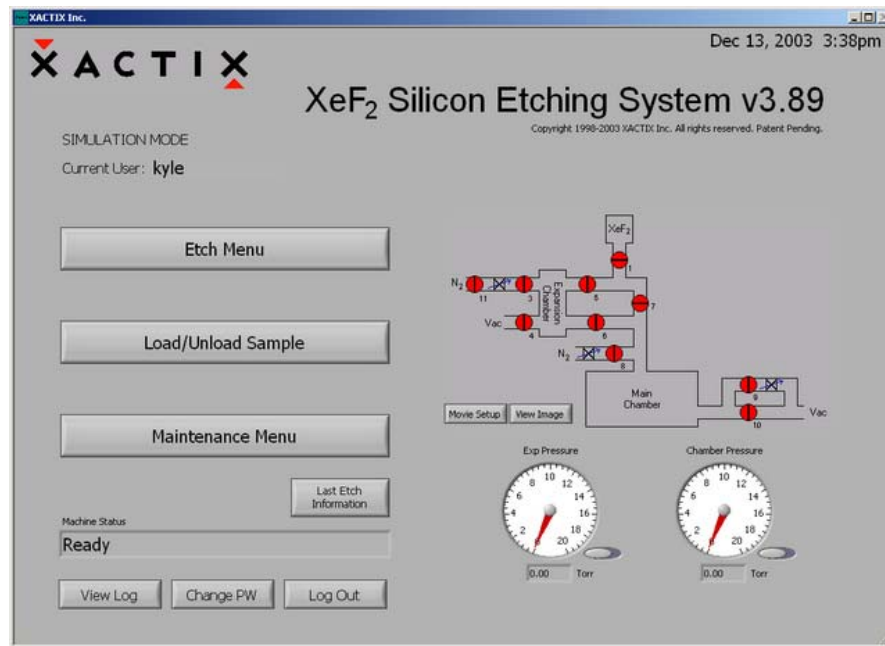


- Press “Load/Unload Sample”. Press “Yes” when prompted, unless you do not wish to load a sample. The system begins chamber purges and flushing cycles to evacuate the chamber.
- If the chamber ventilation shroud is not pulled forward fully, a request to close the shroud will be displayed before venting the chamber.




- Open the chamber lid when the dialog box below appears. Do not press any key prior to opening chamber.

# University of Minnesota Nano Fabrication Center

## Standard Operating Procedure

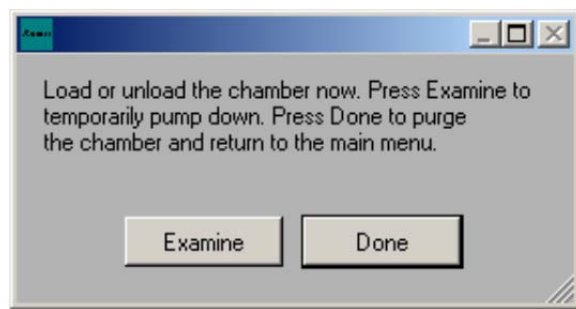

- e. Load a sample. The sample holder and pins should as shown below.

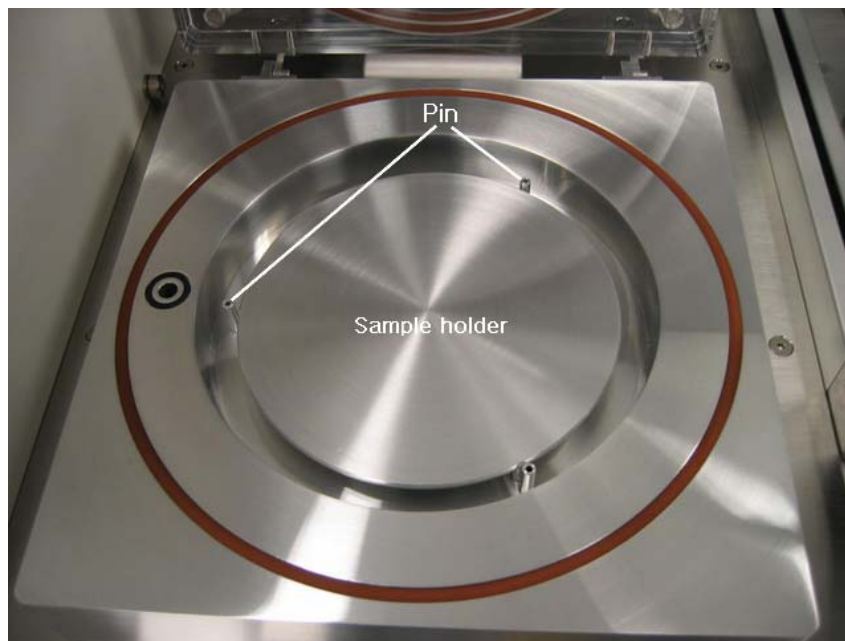


- f. Close lid  
g. Press "Done" on dialog box shown in (c) above. The system will go through a purging cycle prior to chamber pump down.  
h. The Examine button only pumps the chamber down, without purges, so that the system can be quickly vented to load the sample. This is very useful when examining a sample away from the system to prevent moisture from accumulating in the chamber. However, it is always necessary to press DONE before etching the sample.

### Etching

- a. Click on "Etch Menu" in the main menu screen. The screen below appears:  
b. Enter lot number for the sample being etched. You can use alphanumeric characters. Press "DONE".  
c. Two etch operating modes are available:  
1. Normal mode: utilizes a pulsed etch with a set  $\text{XeF}_2$  and  $\text{N}_2$  (if desired) pressure and etch time.  
2. Normal with delay mode: same as normal mode but also allows for a chill time for the sample to cool between each cycle.

# University of Minnesota Nano Fabrication Center

## Standard Operating Procedure

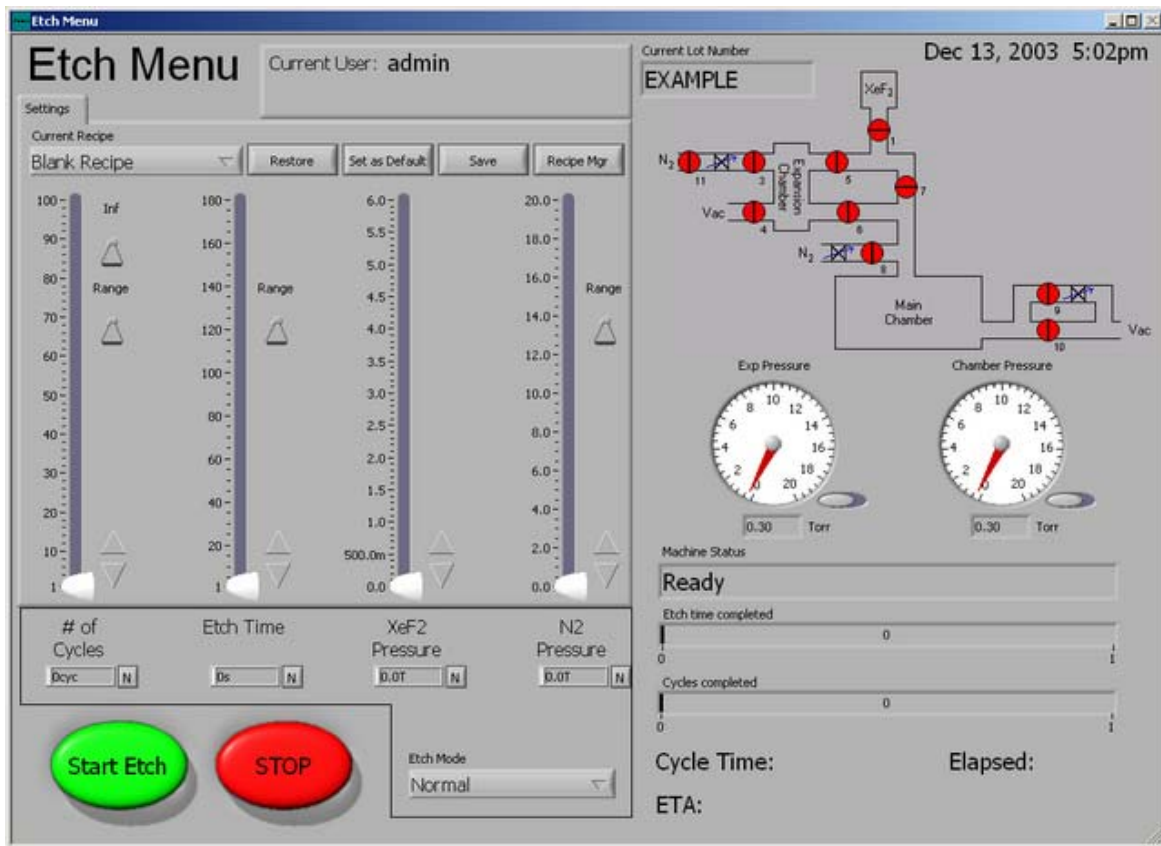

- d. To perform etch in Normal mode: Select normal under “Etch Mode”.
1. Select number of cycles. A cycle consists of the  $\text{XeF}_2$  sublimating to the set pressure in the expansion chamber, etching for a set amount of time, and evacuation of the main and expansion chamber
  2. Select etch time: this time comprises the duration between the opening of the valve between the expansion chamber and the process chamber and the opening of the valve between the process chamber and the pump
  3. Set  $\text{XeF}_2$  pressure: At room temperature, the upper limit for  $\text{XeF}_2$  pressure is 4 torr.
  4. Set  $\text{N}_2$  pressure:  $\text{N}_2$  can be added into a recipe to improve selectivity
- e. Default recipe

# of cycles	Etch time (sec)	$\text{XeF}_2$ (torr)	$\text{N}_2$ (torr)
20	15	1	2


The etch rate for this recipe is about 120nm/min for 50% exposed Si masked with  $\text{SiO}_2$  on a 4-inch wafer.

- f. To perform etch in Normal with delay mode, in addition to the above selections, set the ETCH DELAY which is the time in seconds that the system waits before beginning the next etching cycle.

# University of Minnesota Nano Fabrication Center

## Standard Operating Procedure

- g. Press the “START ETCH” button. Progress of the etch can be monitored following the Etch Time Completed and Cycles Completed bars along with the Elapsed Cycle Time and the ETA.
- h. “CHANGE CYCLES” tab can be used at any time during an etch to adjust the number of cycles. It appears on the screen only when etching has been started. When selected (see dialog box below), use the top row of arrows to add cycles or the bottom row of arrows to remove cycles.

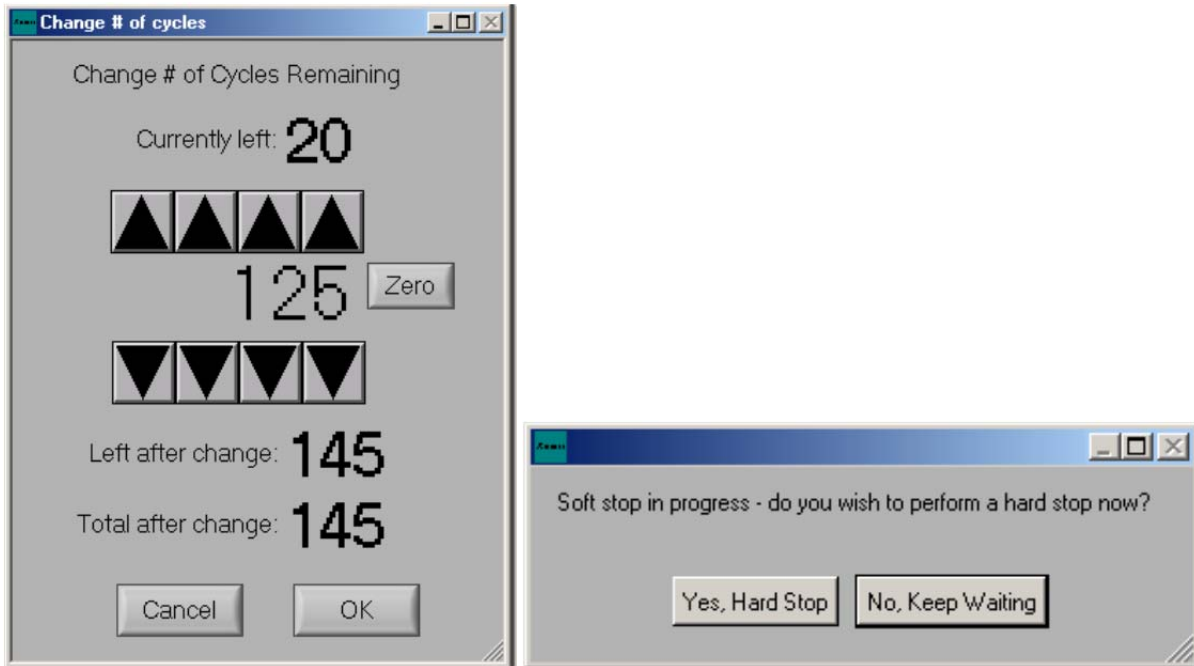

- i. Press STOP to end the etching. The current cycle will be completed before ending the entire etch. If the “STOP” button is pressed twice, the Xetch will prompt the above. Choosing “Yes, Hard Stop” will end the etch cycle immediately and return the user to the main menu. If “No, Keep Waiting” is chosen only the current cycle will be completed before returning the user to the main menu.

### Unloading samples

- a. Press LOAD/UNLOAD and follow the steps as in 8 above.
- b. Unload your samples
- c. Press DONE to pump down the chamber.

### Viewing Detailed Etch Information

- a. The Detailed Etch Information log may be viewed by clicking on the VIEW LOG button on the Main Menu.
- b. The log file is a database that is queried by beginning date and ending date. The “Today” button will automatically set the dates to the current month, day, and year. The lot number, username, recipe, note’s keyword(s) and/or etching mode used can further



# University of Minnesota Nano Fabrication Center

## Standard Operating Procedure

specify your search. Wildcard characters (\*,?) may be used to fully specify the search criteria.

### Creating or Changing a Recipe

- Creating or changing recipes is done from the ETCH MENU screen.
- Make the appropriate recipe changes
- Save the recipe by clicking on the save button located near the top of the “perform etch screen”.
- The save new recipe screen will appear.



- Global Recipes are starting recipe templates available to all users.
- Using the toggle switch in the middle of the screen users can save the new recipe to their Personal Directory or the accessible Global Recipes.
- Type in the name of your new recipe and Click on the SAVE button.